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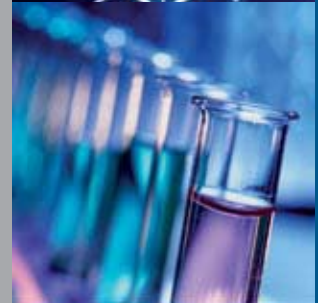
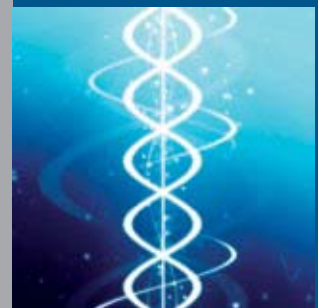
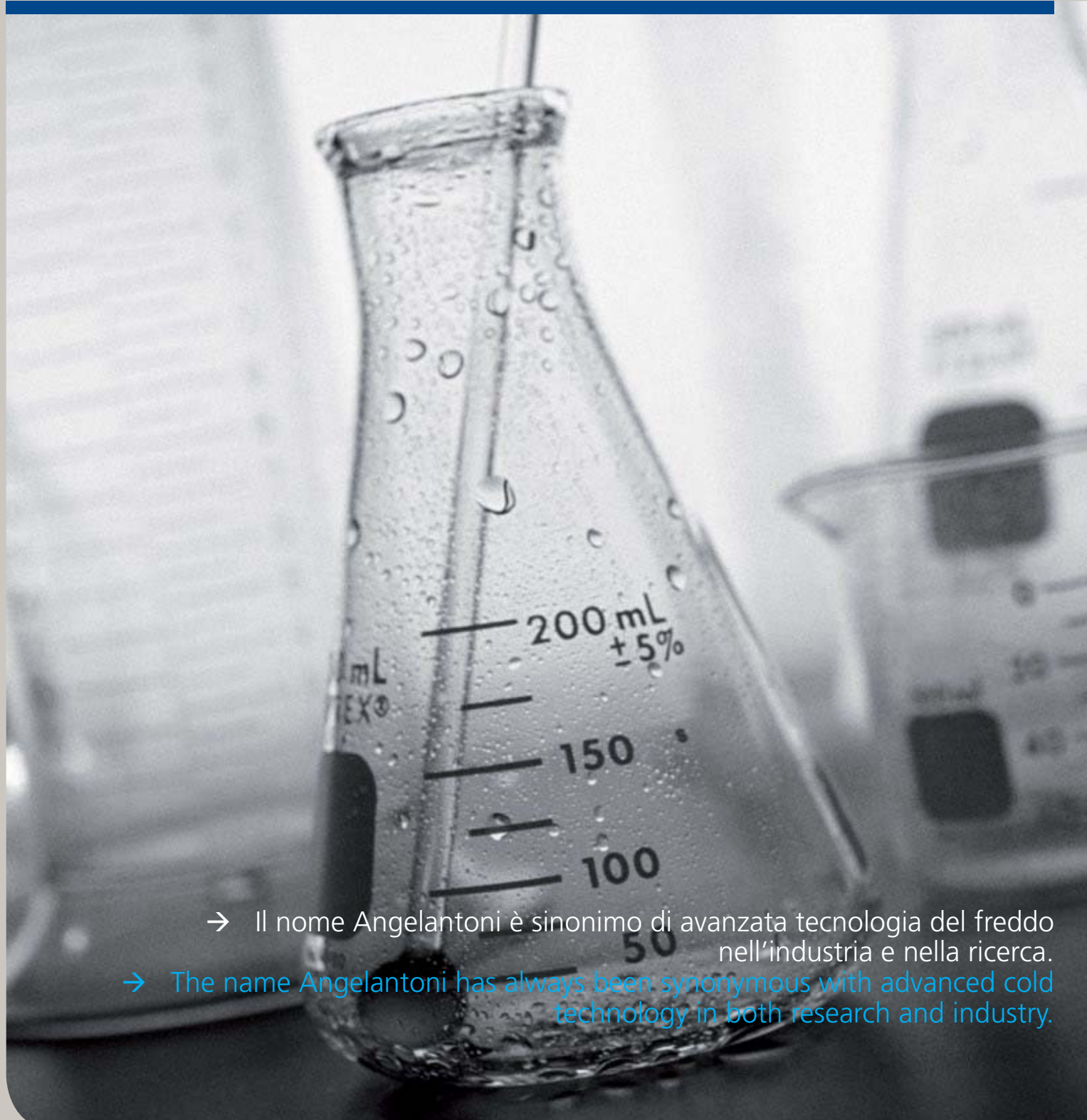
Angelantoni Industrie began operating in 1932 in the refrigeration field, and over the years it has become increasingly specialized in the biomedical sector with the technological and scientific applications of refrigeration, but also in the broader field of laboratory equipment, such as laminar flow hoods and controlled contamination environments.

- **Today Angelantoni Industrie S.p.A. (Holding), comprises three separate subholdings:**

**Angelantoni Life Science S.r.l. (ALS)** with the main brands AHSI – AS – STERIL – AG – ACOTEC (biomedical equipment, controlled contamination environment equipment, laminar flow hoods and clean rooms, instruments and furnishings for research laboratories, health facilities and the pharmaceutical industry);

**Angelantoni Test Technologies S.r.l. (ATT)** with the main brands ACS – TIRA – BIA – AMEC – AKI (environmental test chambers and space simulators, test benches for the automotive industry, electrodynamic shakers, balancing systems);

**Angelantoni CleanTech S.r.l. (ACT)** with the main brands SOLARLIGHT – 3RAYS – KENOSISTEC – ELIANTO – ENTERPRISE (photovoltaic plants for **asbestos replacement, photovoltaic greenhouses, concentration photovoltaic systems, solar thermal power plants with Fresnel reflectors, equipment for thin film technology coating, biogas power plants**).

Associated with ACT, but with a separate structure, ARCHIMEDE SOLAR ENERGY designs and produces molten salt solar receiver tubes, superheated oil and steam (DSG) for solar thermal power plants with parabolic or flat (Fresnel) reflectors.

- **Each of the brands in ALS has many years of experience in the biomedical sector.**

**Angelantoni Scientifica (AS)**. The field of refrigeration equipment for the maintenance and storage of biological materials has received a considerable boost in recent years through Research & Development activities. In 1961 Angelantoni produced the first freezer in Europe able to

reach  $-100^{\circ}\text{C}$  without the use of cryogenic gases. Along with the traditional laboratory refrigerators and the new generation of  $-86^{\circ}\text{C}$  freezers, with the AS brand Angelantoni Life Science has put extremely innovative products on the market, with completely automated control systems, such as the HEMOSAFE™ blood storage systems and the SMARTFREEZER™  $-86^{\circ}\text{C}$  and  $-180^{\circ}\text{C}$  freezers for biological materials complete with **RFID tracking systems**.

**ACOTEC**, an engineering company, is specialized in turnkey setups for controlled contamination and biological containment environments (“clean rooms”), for the pharmaceutical, hospital and microelectronics industries.

**AG**, a historic brand, utilizes its considerable experience in the refrigeration field to design and build refrigeration systems for industrial processes, especially in the chemical-pharmaceutical sectors.

**STERIL**, which is proud of the many national and international awards it has received for its continuous innovation, is the ALS division dedicated to the designing and manufacturing of horizontal and vertical laminar flow cabinets, biohazard and cytostatic safety cabinets, laminar flow pass boxes with UV, sanitizing pass boxes with hydrogen peroxide, cabinets for weighing, sampling and dispensing and isolators for active powders in compliance with the most recent international standards (cGMP).

**AHSI** instead offers to the Italian market a complete range of biomedical and pharmaceutical products and services. Together with ALS equipment, AHSI also offers laboratory furniture, chemical hoods, laboratory centrifuges, CO<sub>2</sub> incubators, nitrogen peroxide sanitizing systems, also available in packages, as laboratory turnkey projects.

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→ Today the ANGELANTONI group includes 6 production units located in Italy, Germany, France, India and China, with a total of over 850 employees and turnover of 145 million Euro.

<div data-bbox="325 443 405 479" data-label="Text">/01. A</div> <div data-bbox="325 786 405 822" data-label="Text">/01. B</div> <div data-bbox="325 1178 405 1214" data-label="Text">/01. C</div> <div data-bbox="325 1294 405 1330" data-label="Text">/01. D</div>	<b>Frigoriferi e congelatori /Refrigerators &amp; freezers</b>	8	
	Temperature range [+4 °C /-30 °C]	<ul style="list-style-type: none"> <li>• UBR/UBF</li> <li>• FRL</li> <li>• FRL PRO</li> <li>• EKOZERO</li> <li>• EKOBASIC</li> <li>• CE</li> <li>• FCL</li> <li>• FCL PRO</li> <li>• EKOFRIGOLAB</li> </ul>	
	<b>Congelatori, ultracongelatori e criocongelatorimeccanici /Freezers, ultrafreezers &amp; mechanical cryofreezers</b>	13	
	Temperature [-30 °C]	<ul style="list-style-type: none"> <li>• KRYOLAB</li> <li>• KRYOLAB PRO</li> </ul>	
Temperature [-45 °C]	<ul style="list-style-type: none"> <li>• KRYOS</li> </ul>		
Temperatures [-40 °C] [-86 °C]	<ul style="list-style-type: none"> <li>• PLATILAB</li> <li>• IRILAB</li> </ul>		
Temperature [-152 °C]	<ul style="list-style-type: none"> <li>• RHENIUM 340</li> </ul>		
<b>Frigoriferi e congelatori portatili /Portable refrigerator &amp; freezers</b>	16		
Temperature range [+10 °C /-18 °C]	<ul style="list-style-type: none"> <li>• MPRF</li> </ul>		
<b>Produttori di ghiaccio /Ice makers</b>	16		
<ul style="list-style-type: none"> <li>• CA    Cubetti /Ice Cubes</li> <li>• FB    Scaglie /Flakes</li> </ul>			
Medical Devices Line	/02. A	<b>Friogemoteche /Blood bank refrigerators</b>	17
<b>93/42/EEC</b>	Temperature [+4 °C]	<ul style="list-style-type: none"> <li>• HEMOSAFE™</li> <li>• BBR</li> <li>• MINI</li> </ul>	
/02. B	<b>Frigoriferi e congelatori prefabbricati /Prefabricated blood bank refrigerators &amp; Freezers</b>	19	
<ul style="list-style-type: none"> <li>• Armadi /Reach-in refrigerators &amp; freezers [+4 °C]</li> <li>• Camere /Walk-in refrigerators &amp; freezers [-30 °C]</li> </ul>	Orizzontali e verticali / Horizontal & vertical [-45 °C] [-86 °C]		
<ul style="list-style-type: none"> <li>• KRYOS</li> <li>• PLATINUM™</li> <li>• IRIDIUM™</li> </ul>			
/02. C	<b>Congelatori rapidi per plasma /Plasma blast freezers</b>	22	
Temperature [-70 °C]	<ul style="list-style-type: none"> <li>• PLASMAFROST™</li> <li>• PLASMAFROST NEXT ITEM™</li> </ul>		
/03. A	<b>Frigoriferi e congelatori per latte materno /Mother milk refrigerators &amp; freezers</b>	24	
<ul style="list-style-type: none"> <li>• MMR [+4 °C]</li> <li>• MMF [-20 °C]</li> </ul>			

/04. A	<b>Camere climatiche, termostatiche e fitologiche /Stability test and plant growth chambers</b> 25 <ul style="list-style-type: none"><li>• EKOCT EKOCTL [+10 °C /+50 °C]</li><li>• EKOCH EKOCHL [-10 °C /+60 °C]</li></ul>
/05. A	<b>Anatomia patologica - Medicina legale /Morgue equipment and mortuary chambers</b> 27 <ul style="list-style-type: none"><li>• PM TN (Medium conservation) [+4 °C]</li><li>• PM BT (Long conservation) [-20 °C]</li><li>• Carrelli e accessori /Accessories and carts</li><li>• EG</li></ul>
/06. A	<b>Soluzioni software per laboratorio /Laboratory software solutions</b> 29 <ul style="list-style-type: none"><li>• FREEZEPRO</li><li>• SPYLOG PLUS</li></ul>

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Opening system based on a transponder key, opening door occur with a personal magnetic TAG.

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Software to monitoring the equipment through a PC, with the following specs: BIOGUARD® door opening management system; temperature control; microprocessor control board.

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Angelantoni Industrie SpA offers a wide range of refrigerated cabinets at single or dual temperature (combination of a refrigerator and a freezer in a single unit) dedicated to the preservation of biological materials and pharmaceuticals.

In particular, the broad range of models of different capacities fully copes with diverse needs of the healthcare and biotechnology.

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#### 120 and 170 liters under bench reach-in refrigerators:

- UBR series\_ [ +2°C / +12°C ]
- UBF series\_ [ -15°C / -30°C ]

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
Abrasion proof special white plastic internal execution.  
Safety key lock.  
Temperature digital display.  
Mechanical thermostat for temperature regulation.  
N° 4 adjustable shelves, one in glass (UBR only).  
N° 2 aluminum fixed shelves and n° 3 drawers (UBF only).  
Internal lamp (UBR only).

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### Refrigerators from 160 to 500 liters.

- FRL series\_ [ +2°C / +12°C ]

Fully metallic in phosphate-coated, powder-painted steel plate.  
Internal chamber with abrasion proof special white plastic.  
Service door with variable hinges according to needs.  
Safety key lock.  
Grid shelves, adjustable in height.  
Temperature digital display and internal ventilation.

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### Laboratory refrigerators, FRL-PRO series. Capacities: 100, 200 e 300 Liters.

Temperature range: [+2°C /+12°C]

Main structure in galvanized steel anti-corrosive and non-toxic, coated with PVC film gray. Solid door or, on request, double glazed (GL version), self locking, magnetic seal.  
Rollers for easy movement.

Thermal insulation, thickness 60 mm, obtained by the injection of polyurethane foams, HCFC free, under strong pressure ("sandwich technique").  
Inner chamber, equipped with automatic lighting, stainless steel AISI 304 with Scotch Brite shelves. Internal evaporator fan. Hole Ø 23 mm with a rubber stopper.

#### REGULATION AND CONTROL SYSTEM

- Digital-Display.
- Temperature-control microprocessor.
- Turning on and off-device via key switch ON / OFF.

#### AUDIBLE AND VISUAL ALARMS

- High and low temperature alarm.
- Open door alarm.
- Clogged condenser alarm.
- Alarm faulty probe.
- Voltage failure alarm, self-powered control unit to ensure within 48 hours after the operation of the control system.

Possibility of remote alarms. Storing alarms. The system is equipped with two analog inputs for sensors, temperature sensing probe PT1000; KTY sensor for sensing temperature condenser.

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Optional

#### Laboratory refrigerators, EKOZERO series. Available capacities: 700 and 1500 liters.

- **TN series\_** [ 0°C / +15°C ]

Fully metallic powder-painted in heavy-gauge phosphate-coated. AISI 304 Stainless Steel internal executions with rounded corners. Average thickness of the insulating layer: 75 mm. Safety key lock. Reversible door, made of the same features of the main structure, with special magnetic seals. Reversible door in triple glazed is available on request. Automatically locking system for door open beyond 90°. Height adjustable feet. Internal light, with manual or automatic switch. Internal ventilation. Height adjustable shelves. Maximum external temperature +32°C.

Electronic control with:

- Digital display indicating the temperature inside the refrigerator cabinet.
- Power on / off camera.
- Key to activate manual defrost.
- Button setting the temperature.
- Manual activation button for internal light.
- Visual and audible alarm for open door, low and high temperature.
- Storage of minimum and maximum temperature reached.
- Preparing the remote alarm output, with NA 230V contact.

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Optional

#### 700 and 1500 liters Laboratory Refrigerators:

- **TN series\_** [ 0°C / +15°C ]
- **BT series\_** [ -10°C / -20°C ]

Fully metallic powder-painted in heavy-gauge phosphate-coated. AISI 304 Stainless Steel internal executions. Average thickness of the insulating layer: 75 mm. Safety key lock. Port hole 23 mm. Internal ventilation. Internal light, with manual or automatic switch. Grid shelves, adjustable in height. Microprocessor electronic control system. Remote alarms.

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Contatti allarme remoto.

**Modular prefabricated Refrigerators with capacity from 2000 to 5000 liters:**

- **TN series\_** [ 0°C / +10°C ]
- **BT series\_** [ -10°C / -20°C ]

Walls, floor and ceiling in prefabricated modular panels:

- Thickness of panel: 60 mm.
- External and internal layer in white prepainted zinc plate steel.

Safety key lock.

Internal ventilation.

Internal light, with manual or automatic switch.

Grid shelves, adjustable in height.

Proprietary microprocessor control system (using PT100 probe).

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N° 2 ruote posteriori e n° 2 piedini anteriori.

Chiusura con chiave esterna.

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**Refrigerators /Freezers with two independent chambers and cooling units of 300 and 400 liters.**

- **FCL series\_** [ Refrigerator: +1°C /+9°C.  
Freezer: -15°C / -32°C ]

Fully metallic in phosphate-coated, powder-painted steel plate.

Internal chamber with abrasion proof special white plastic.

Average thickness of the insulating layer for the refrigerator chamber: 40 mm.

Average thickness of the insulating layer for the freezer chamber: 80 mm.

N° 2 back rollers and n° 2 front rollers.

External lock with key.

Internal light (+4°C compartment only).

Internal ventilation (+4°C compartment only).

Shelves and drawers in the refrigerator chamber.

Drawers in the freezer chamber.

Temperature led indicator for each compartment.

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**Refrigerator / freezers with two independent compartments with dual refrigeration system, from 300 to 700 liters.**

**FCL-PRO Series**

**Refrigerator:** [+2°C /+12°C]

**Freezer:** [-10°C /-25°C]

Main structure in galvanized steel anti-corrosive and non-toxic, coated with PVC film gray.

Insulation thickness: 60 mm.

Rollers for easy movement.

N° 2 back doors self-closing metal painted gray, complete with special magnetic strips, electrical resistance heater placed on the frames of doors, with lock and key. On request the door of the refrigerator compartment is also available in glass.

N° 2 holes (ø 23 mm with a rubber stopper).

Interior lighting (only for +4 °C compartment).

Ventilation system (only for +4 °C compartment).

Compartment at +4 °C:

N° 3 stainless steel shelves adjustable in height.

Compartment at -25 °C:

Drawers and removable stainless steel mounted on telescopic rails.

Optional

### Ekofrigolab series\_ Professional line

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Montaggio su sfere (700-850-1500).

Piedini regolabili (130-250).

**Professional lab refrigerators with capacity from 130, 250, 700, 850 and 1500 liters:**

- **TN series\_** [ 0°C / +15°C ]
- **BT series\_** [ -10°C / -23°C ]
- **LT series\_** [ -10°C / -30°C ]
- **2TS series\_** [ Refrigerator: 0°C /+15°C. Freezer: -10°C / -25°C ]

Fully metallic powder-painted in heavy-gauge phosphate-coated.

AISI 304 Stainless Steel internal executions.

Average thickness of the insulating layer: 75 mm.

Porthole 23 mm.

Safety key lock.

Internal light, with manual or automatic switch

Internal ventilation.

Shelves, adjustable in height.

Proprietary microprocessor control system (using PT100 probe).

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Angelantoni Industrie SpA offers a broad range of freezers at -20°C, -30°C, -45°C and -85°C available to meet the expectations of healthcare and biotechnology. Particularly suitable for use in hospitals and laboratories for freezing and long storage of microorganisms, viruses, or any kind of biological sample. IRILAB<sub>Next</sub> represents the new line of freezers at -80 /-85°C, able to develop functionality, performance and correct preservation of any biological material.

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**Chest and upright freezers with capacity from 300 to 500 liters.**

• **Kryolab series\_** [-15°C / -32°C ]

Fully metallic in phosphate-coated, powder-painted steel plate.  
 Internal chamber with abrasion proof special white plastic.  
 Safety key lock.  
 Temperature LED indicator for KRYOLAB 300H, 500H and 300V.  
 Digital display temperature indicator-regulator KRYOLAB 500V.  
 Drawers and shelves.

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**Upright freezers KRYOLAB 320 V-PRO Series.**

**Temperature range:** [-10°C / -30°C ]

External structure of sheet zinc hot plastic anti-corrosion and non-toxic in the color gray.  
 N° 1 service port of the same characteristics of the structure with stainless steel interior, complete with special magnetic strips, electrical resistance heater placed on the door frame, lock and key. Rollers for easy movement.  
 Porthole Ø 23 mm with a rubber stopper.  
 Internal chamber in stainless steel AISI 304 with rounded corners.  
 Insulation, 80 mm, obtained by the injection of polyurethane foams, HCFC free, under strong pressure ("sandwich technique").  
 N° 3 sliding drawers and removable stainless steel, mounted on telescopic guides which allow the complete discharge of the drawer.  
 Regulation system by digital display and control by microprocessor.

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**Horizontal and vertical freezers of 500 liters. Series Kryos.**

- Kryos SH series [ -25°C / -45°C ]
- Kryos SV series [ -20°C / -43°C ]

Structure in sheet steel epoxy coated.

Internal chamber in aluminum (SH model) or in white special anti-abrasion plastic (SV models).

Internal lighting (SH model).

Door with key (STD version), RFID system with transponder (Bioguard® PLUS version, available only on SV models).

Microprocessor digital temperature regulator.

Remote alarm.

Baskets (SH model), n. ° 14 drawers (SV model).

Optional

### Platilab™ series\_ Professional line

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Montaggio su ruote.

**Chest and Up-right freezer line from 113 to 550 liters.**

- SH-SV series [ -45°C ]
- H-V series [ -86°C ]

Fully metallic in heavy-gauge phosphate-coated, powder-painted steel plate.

Safety key lock.

AISI 304 Stainless Steel internal execution.

Insulated inner doors.

Stainless steel internal shelves.

23 mm Port hole.

Proprietary microprocessor control system (using PT100 probe).

Mounted on wheels.

Optional

## Irilab<sub>Next</sub> series\_ Professional line

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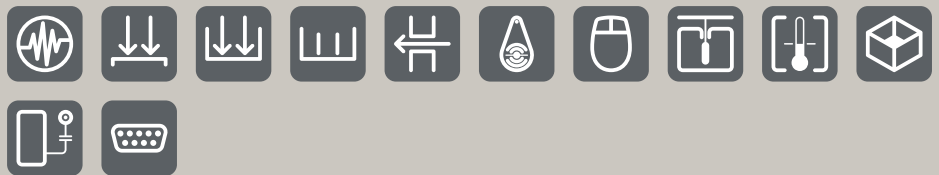
### 520 and 800 liters up-right freezer.

- V series\_ [ -80 /-86°C ]
- SV series\_ [ -45°C ]

10/10 AISI 304 Stainless Steel internal execution.  
Mounted on wheels.  
NOT insulated inner doors.  
Magnetic lock with key.  
Stainless steel internal shelves.  
23 mm Port hole.  
New Microprocessor control system COLD BRAIN with graphic display 240x128 pixels, which implements in a single video "Regulation and Alarms Section" using 8 soft keys.  
Ethernet interface (Wireless Wi Fi optional).

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### Mechanical Cryofreezer at -152 ° C with a capacity of 340 liters.

- Rhenium series\_ [ -135°C /-155°C ]

#### RHENIUM: THE LARGER STORAGE CAPACITY MECHANICAL CRYOGENIC FREEZER IN THE MARKET!

External plate frame, painted with grey epoxy resins. N° 1 phosphate steel plate door, painted with grey epoxy resins, with magnetic gaskets and heavy duty handle. The door is equipped with key lock (STD version) or RFID transponder system with Bioguard® (PLUS version).

N° 6 wheels to facilitate movement of the freezer.  
Internal chamber in stainless steel 8/10 AISI 304; rounded corners for easy cleaning.  
Thermal insulation with high thermal efficiency obtained with polyurethane foam and vacuum insulation panels (VIP).  
Insulation thickness: 140 mm.

New control system with COLD BRAIN user interface with graphical display 240 x 128 pixels, which implements a single video on the "Regulator Section" and "Box Alarm" through a capacitive touch keypad (8 soft keys).

Serial interface standard (wireless option).

Optional

Angelantoni Life Science /Range of production AS /01. C-D

/01. C **Frigoriferi e congelatori portatili** /Portable refrigerator & freezers

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Angelantoni Industrie SpA portable refrigerators provide a safe transportation of biological samples and blood products and ensure the best storing conditions.

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#### **Portable refrigerators and freezers.**

Electronic thermostat temperature regulator.  
Visual temperature display.  
AC /DC current.  
Double voltage: **DC** 12 /24 and **AC** 115 /230 V.

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Angelantoni Industrie SpA manufactures ice cubes and flakes makers of high quality to guarantee the best performance.

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#### **Complete range of ice maker or cube makers with daily variable output.**

Shockproof plastic doors.  
A safety thermostat is installed on the equipment to stop it automatically once the maximum level of ice in the container is reached.

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Angelantoni Industrie SpA has a wide range of blood bank refrigerators (BBR) certified according to Directive 93/42/EEC and all modifications, to meet demands of small, medium and large transfusion centers for hospitals, analytical laboratories and operating rooms. The models available can store 30 to 1.400 blood bags and ensure, through a proprietary electronic system, the internal temperature at +4°C, as required by the current international standards.

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#### Automatic blood bank refrigerators/ Hemosafe®

HEMOSAFE® is the most safe blood bank refrigerator in the world: the right blood to the right patient at the right time!

HEMOSAFE® is the only automated and computerized blood bank refrigerator for a safe and smart management of blood units stock (up to 150).

Thanks to the integrated management software, HEMOSAFE® is able to select the most suitable blood bag for each patient and to deliver it to the operator quickly (in less than 60 seconds!) and without any human mistakes. Operator access is controlled and limited to the right blood bag.

HEMOSAFE® makes the blood available 24/7 in each remote hospital, optimizing the distribution and reducing the number of journeys between Blood Center and hospitals.

HEMOSAFE®, with Just in Time delivery and Electronic Crossmatch, guarantees a significant reduction in the number of unused blood units and eliminates unnecessary labeling and pre-allocation activities. FIFO (First In First Out) logic allows an ideal stock management according to expiration dates.

HEMOSAFE® can be easily interfaced with any LIMS (Laboratory Information Management System) and BBIS (Blood Bank Information System).

HEMOSAFE® follows the best practices and recommendations by the most eminent transfusion associations in the world (AABB, ISBT, SIMTI) about the implementation of Type and Screen and Electronic Crossmatch.

HEMOSAFE® is a Medical Device certified according to 93/42/EEC and following amendments.

**Capacity:** 150 blood bags (450)  
**Temperature:** +4°C  
**Size (External) mm:** 1045 x 890 x 1845 (W,D,H)

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#### Monolithic or prefabricated blood bank refrigerators.

- BBR series\_ [ from 200 to 1064 bags ]

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
 Full-view glass doors.  
 AISI 304 Stainless steel internal execution.  
 Safety key lock.  
 Average thickness of the insulating layer: 75 mm.  
 Port hole 23 mm.  
 Internal light, with manual or automatic switch.  
 Internal ventilation.  
 Proprietary microprocessor control system (using PT100 probe).  
 Remote alarm contacts.  
 Tropical execution.  
 Temperature graph recorder.  
 Stainless steel sliding drawers.  
 Mounted on spherical wheels.

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#### Monolithic blood bank refrigerators.

- Mini series\_ [ from 30 to 100 bags ]

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
 Full-view glass doors.  
 Zinc plated steel pre-painted in white internal execution.  
 Safety key lock.  
 Average thickness of the insulating layer: 40 mm.  
 Port hole 23 mm.  
 Internal light, with manual or automatic switch.  
 Internal ventilation.  
 Proprietary microprocessor control system (using PT100 probe).  
 Remote alarm contacts.  
 Temperature graph recorder.  
 Stainless steel sliding drawers.

Optional

Notified Body No.:  
0434

## /02.B

### Armadi e camere / Reach-in & Walk-in refrigerators & freezers

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#### Prefabricated blood bank refrigerators and freezers PE Series.

Reach-in refrigerators and freezers (from 3000 to 6000 liters).

- **TN models** \_ [0°C / +10°C]
- **LT models** \_ [-10°C / -30°C]

Zinc steel plate main frame made of white pre-painted prefabricated panels with gray upper strip. Up to 6 service doors with handles, locks and specific rubber gaskets. Anti-mist electric resistances on the doors frame. External key lock. Inner chamber in abrasion proof special white plastic. Evaporator placed on fixed shelves. Up to 24 shelves.

Isolation obtained by "sandwich technique", with polyurethane foams. Insulation gauge: 60 mm (TN models); 120 mm (LT models).

Adjustment and control system by E2003 microprocessor, with two independent sections, the first one to set and control temperature, the second one to manage the alarm system and display temperature.

#### Walk-in chambers (from 10000 to 21000 liters).

- **TN models** \_ [0°C / +10°C]
- **LT models** \_ [-10°C / -30°C]

FLOOR / Stainless steel coated with anti-glaciation grids. Prefabricated chamber made of modular self-holding welding less panels in order to allow a total disassembly.

The door is in the front wall.

Adjustment and control system by microprocessor controller E2003, with two independent sections, the first one to set and control temperature, the second one to manage the alarm system and display temperature.

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Complete range of freezers at -40°C certified according to Directive 93/42/EEC and all modifications. These devices are specifically adapted for use in hospitals and laboratories for freeze and storage in long periods, blood components and derivatives.

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#### Vertical freezers of 500 liters. Series Kryos.

- **Kyos SV Series** [-20 °C / -43 °C]

Structure in sheet steel epoxy coated.  
Internal chamber in white special anti-abrasion plastic with n. ° 14 drawers.  
Door with key (STD version), RFID system with transponder (Bioguard® PLUS version).  
Microprocessor digital temperature regulator.  
Remote alarm.  
Equipment certified according to 93/42/EEC directive and following amendment.

Optional

Platinum™ series\_ Professional line

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Chest and Up-right freezer line from 113 to 550 liters.

- SH-SV series\_ [ -45°C ]
- H-V series\_ [ -85°C ]

Fully metallic in heavy-gauge phosphate-coated, powder-painted steel plate.  
Safety key lock.  
AISI 304 Stainless Steel internal execution.  
Insulated inner doors.  
Stainless steel internal shelves.  
23 mm Port hole.  
Proprietary microprocessor control system (using PT100 probe).  
Mounted on wheels.

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800 liters up-right freezer.

- SV series\_ [ -40°C ]
- V series\_ [ -80°C ]

Rigid box type frame, powder-painted steel plate.  
AISI 304 Stainless Steel internal execution.  
NOT insulated inner doors.  
Magnetic lock with key.  
Stainless steel internal shelves.  
23 mm Port hole.  
Proprietary microprocessor control system (using PT100 probe).  
Mounted on wheels.

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Specifically built to freeze the blood plasma at  $-30^{\circ}\text{C}$  in less than 60 minutes in compliance with the most restrictive standards to preserve the plasma active ingredients and certified according to Directive 93/42/EEC and all modifications.

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#### Blast Freezers for plasma bags.

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
AISI 304 Stainless steel internal execution.  
Service door with key lock.  
Average thickness of the insulating layer: 125 mm.  
Shelves and uniformity plates.  
Mounted on wheels.

Control panel:

- On /Off button.
- Signal light for cooling phase.
- Visual and acoustic signal for cooling cycle.
- Shelf temperature digital display.
- On/Off main switch.

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### Plasmafrost ITEM™ series.

The plasma blast freezers Plasmafrost I.Te.M. series (Indirect Temperature Measurement) monitor each cycle of freezing processes and certify that the inner part of each plasma bag reach a temperature of -30 °C in less than 60 minutes, as required by the European recommendation R (95) 15.

The core of the technology is the system I.Te.M., that by measuring the temperature of the plate of uniformity and through the use of an algorithm implemented in software in the integrated touch-screen PC, makes possible the reliable and precise temperature detection of the interior of each bag of plasma identified with the bar codes reader (integrated) and loaded into the freezer.

The report that Plasmafrost I.Te.M. print at the end of the freezing process include the following details:

- Starting and final freezing times
- Plasma bags barodes
- Process steps details
- Operator ID
- Report date and time
- Signature of the validating officer

Plasmafrost I.Te.M. thanks to the touch-screen panel PC, barcode reader and intgrated proprietary software allows:

- To identify of all the bags in a freezing batch
- To measure the temperature of each bag in a freezing batch
- To record and back trace the freezing process data
- To protect the freezing data from manipulation
- To share the freezing data with hospital/blood center information systems

Angelantoni Life Science /Range of production AS /03. A  
Frigoriferi e congelatori per latte materno  
Mother milk refrigerators & freezers

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The MMR and MMF refrigerators and freezers from 700 to 1500 liters are ideal to store human milk, a key product for the development and future health of the children. These equipment are managed by dedicated software to secure the correct distribution of milk and its proper administration to premature babies in intensive care.

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**700 and 1500 liters Mother Milk Refrigerators and Mother Milk Freezer** to perfectly preserve and store the mother milk.

- MMR series\_ [ 0°C / +15°C ]
- MMF series\_ [ -10°C / -23°C ]

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
Full-view glass doors.  
AISI 304 Stainless steel internal execution.  
Bioguard® access control system.  
Average thickness of the insulating layer: 75 mm.  
Port hole 23 mm.  
Internal light, with manual or automatic switch.  
Internal ventilation.  
Proprietary microprocessor control system (using PT100 probe).  
Remote alarm contacts.  
Tropical execution.  
Temperature graph recorder.  
Mounted on spherical wheels.

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Stability tests at controlled temperature and relative humidity according to current GLP and GMP standards are gaining increasing importance in drugs and cosmetics stability. Stability tests at controlled temperature, relative humidity and light are used in biotechnology, entomology, botany and the food industry. The rather small external dimensions, easiness of use, reduced energy consumption and high reliability are the success keys of this line.

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**700 and 1.500 liters thermostatic chamber.**  
 Range of temperature: from +10°C to +50°C.

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
 Stainless steel internal execution.  
 Glass inner door.  
 Safety key lock.  
 Average thickness of the insulating layer: 75 mm.  
 23 mm port hole.  
 Internal ventilation.  
 Stainless steel AISI 304 grid shelves with supports, adjustable in height.  
 Temperature electronic regulator at digital reading.  
 Mounted on spherical wheels.

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**700 and 1.500 liters solar simulation chamber** enable the combination of temperature cycles with indoor lighting. They provide variable lighting of 20.000 lux.

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
 Stainless steel internal execution.  
 Glass inner door.  
 Safety key lock.  
 Average thickness of the insulating layer: 75 mm.  
 23 mm port hole.  
 Internal ventilation.  
 Stainless steel AISI 304 grid shelves with supports, adjustable in height.  
 Temperature electronic regulator at digital reading.  
 Relative Humidity electronic regulator.  
 Lighting system control is carry out by 7 - days timer and On-Off switch (light intensity 0 or 20.000 lux).  
 Mounted on spherical wheels.

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#### Complete range of climatic chamber of 700 and 1.500 litres for drugs stability testing.

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
 Stainless steel internal execution.  
 Glass inner door.  
 Safety key lock.  
 Average thickness of the insulating layer: 75 mm.  
 23 mm port hole.  
 Internal ventilation.  
 Stainless steel AISI 304 grid shelves with supports, adjustable in height.  
 Temperature electronic regulator at digital reading.  
 Relative humidity microprocessor control regulator at digital reading.  
 Mounted on spherical wheels.

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#### 700 and 1.500 liters solar simulation chamber enable the combination of temperature and humidity cycles with indoor lighting. They provide variable lighting of 20.000 lux.

Fully metallic powder-painted in heavy-gauge phosphate-coated.  
 Stainless steel internal execution.  
 Glass inner door.  
 Safety key lock.  
 Average thickness of the insulating layer: 75 mm.  
 23 mm port hole.  
 Internal ventilation.  
 Stainless steel AISI 304 grid shelves with supports, adjustable in height.  
 Temperature electronic regulator at digital reading.  
 Relative humidity microprocessor control regulator at digital reading.  
 Lighting system control is carry out by 7 -days timer and On /Off switch (light intensity 0 or 20.000 lux).  
 Mounted on spherical wheels.

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A quite broad range of products, from mortuary prefabricated rooms of different capacities and temperature (as needed for the conservation of corps) to mortuary complementary articles which allow the transport of the corpses and autopsy tables available in different versions: fixed, elevating and rotating.

### Camere mortuarie /Mortuary chambers PM series\_

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**Standard prefabricated mortuary chamber; medium and long term preservation**, in special execution with body lifting.  
 Turnkey units for morgues.

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- Thickness of panel: 80 mm.
- External and internal layer in white pre-painted zinc plate steel.

Single door with key lock.  
 Stainless steel internal floor.  
 Stainless steel stretchers.  
 Internal light.  
 Telescopic sliders for stretchers.  
 Proprietary microprocessor control system (using PT100 probe).

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**Stainless steel autopsy tables: fixed, stationary and rotating models.**

18/10 Stainless steel.  
 Rounded corners for easy cleaning, disinfection and safety.  
 Support perforated grid for table top.  
 Movable shower fixed to the central basement to supply warm and cold water to the working surface.  
 Table washing system by means of nozzles mounted on the circumference of the table top, manually controlled (not for EG 104).  
 Suction system for liquids (not for EG 104).  
 Suction system for fumes and table air recycling with centrifugal (not for EG 104).

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#### Dissection tables.

18/10 Stainless steel.  
Rounded corners for easy cleaning, disinfection and safety.  
Support perforated grid for table top.  
Movable shower fixed to the central basement to supply warm and cold water to the working surface.

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#### Optional and carts to move and display bodies.

##### Carrier transporting one stretcher, model CTB 1.

Supporting structure entirely made of 18/8 AISI 304 stainless steel.  
Surface support of the stretcher with guides.  
Shackles for the carrier handling included.  
Four pivoting wheels to facilitate movement and blocking clamps.

##### Carrier transporting two stretchers, model CTB 2

Supporting structure entirely made of 18/8 AISI 304 stainless steel.  
Surface support of the two stretchers with guides.  
Shackles for carrier handling included.  
Four pivoting wheels to facilitate movement and blocking clamps.

##### Corpse exhibition table, model TES 1.

Supporting structure entirely made of 18/8 AISI 304 stainless steel.  
Inclined surface with perimetric edges.  
Sewage disposal hole with a removable collection container.

##### Corpse exhibition table, model TES 2.

Made with the same characteristics as the above-mentioned model, but equipped with four pivoting wheels with blocking clamps.

##### Carriers for transport & hoisting of stretchers, model CTS 1.

Supporting structure made of coated steel. Hoist able to the required height by means of a hydraulic piston driven by a suitable pump and easily operated by pedal.  
Equipped with sliding rollers on bearings for easy loading of the stretcher.  
The first roller is equipped with a stopper to prevent movement of the stretcher during transportation.  
Four pivoting wheels to facilitate the movement, the anterior ones with blocking clamps.

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Frozen sample management solution FreezePro® series.

- FreezePro® Standard Version: entry level
- FreezePro® Enterprise Version: business level

FreezePro 2012 is a class-leading frozen sample management solution that is indispensable to any modern scientific or pharmaceutical laboratory.

FreezePro is an intuitive, fast, reliable and secure Web-based application that has a visually appealing interface which enables users to know precisely where a frozen sample is located even before opening the freezer door. Popular features include automatic alerts of low number of sample aliquots, sample expiration date, sample volume or freeze-thaw count along with powerful reporting. FreezePro could be interfaced with RFID or bar code labels for the automatic management of the samples.

Independently validated and certified to be fully compliant with the FDAs current GLP/GMP requirements FreezePro takes laboratory record management to the highest level of productivity!

FreezePro is compliant with the following standards:

- FDA 21 CFR Part 11.
- cGLP/cGMP (Good Laboratory Practice and Good Manufacturing Practice).
- Section 508 Accessibility Guidelines.
- HIPAA Accountability Act of 1996.
- HIPAA Patient Safety and Quality Improvement Act of 2005 (Patient Safety Act).

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### Wireless monitoring system Spylog PLUS

SpyLOG PLUS will keep constantly monitored refrigerators, freezers, blood banks, chambers and environments, where the precious samples and drugs are stored, as well as cold rooms, incubators and other equipment in which the chemical/physical parameters are key for obtaining results.

Several chemical-physical parameters can be acquired and monitored:

- Temperature.
- Relative humidity %.
- Pressure.
- Environments and instruments values of O2%, CO2%, N%.
- Air flow velocity.
- Measures of absolute and differential pressure.
- Environment measures of volatile gas.
- Lighting value (Lux).

#### FIELD OF APPLICATION

The usual appliance is the control of environments and equipment in the following sectors:

#### Hospitals or clinical institutes:

- Transfusion centers.
- Laboratories of analysis, Microbiology, Pathological anatomy.
- Environmental control of Steril Chambers, Operating-Theatres.

#### Pharmaceutical, Chemical, Food and Aerospace companies:

- Quality Control Laboratories.
- Control of Manufacturing area, Warehouses and Pounds.

#### Universities, Research centers, States of territorial control, Animals Health Institutes, museums:

- Development and research laboratories.
- Environments control.

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Companies interested in pursuing success and growth must focus not only on innovation and technology, but also on environment, safety, and human resource aspects. While ISO 9001 certification guarantees compliance with standards, efficiency of processes, and customer satisfaction, Angelantoni Industrie has introduced the **Integrated Management System**, demonstrating that it is possible to apply the principles and rules of quality, environmental respect, and worker and product safety in the manufacture of its products, focusing constantly on accident prevention, process and product improvement, and customer satisfaction.

Angelantoni Industrie has attained:

- **NATO AQAP 104** in 1991
  - **ISO 9001** in 1995
  - **ISO 14001** in 2001
  - **EMAS** (registration number IT-001058) in 2008
  - Product certificates for HEMONINE 2, HEMOSAFE, IRIDIUM starting from 2008
- We also guarantee:
- from 2001: **SA8000, OHSAS 18001** and **SGLS** standards
  - compliance with Directive **2002/95/EC** (RoHS Directive)
  - **RAEE** registration (registration number IT 08020000003520)

In 2009 Angelantoni Industrie also obtained **ISO 13485:2003** certification and the **CE 93/42** mark for medical devices, with their registration with the Ministry of Health: these certifications require the maintaining of the configuration status of medical device products, updating of the personnel, constant monitoring of the performance of the medical devices, a rigorous analysis of complaints and nonconformities, and a more collaborative relationship with customers for any problems.

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Angelantoni Industrie headquarters in Massa Martana (Perugia, Italy) extend over an area of 80.000 square metres (more than 16.000 covered square meters).

Massa Martana is located in Umbria, a region rich in art, history and tradition.

No location could be more appropriate; Angelantoni Industrie S.p.A. learns from the past to better understand and anticipate the future.

This, combined with dedication and over growing expertise, is why Angelantoni Industrie S.p.A. has become the most complete and diversified European Group for advanced cold technology and test equipment for industry and research.

**Our core competencies and services for total customer satisfaction:**

- Training, both at our facility and at customer site.
- Testing and quality checks.
- Installation and start up.
- Service.
- Calibration using SIT certified instruments.
- "Full risk" assistance contracts.
- Special applications /Turn-key project.

ANGELANTONI  
GROUP  
MORE THAN  
YOU THINK

---

**Angelantoni Life Science S.r.l.**

Località Cimacolle, 464 - 06056 Massa Martana (PG) Italy  
ph. (+39) 075.89551 (a.r.) - fax (+39) 075.8955312  
biomedical@angelantoni.it

**[www.angelantonilifescience.it](http://www.angelantonilifescience.it)**

---